

Start Up LSM 800

1. Turn on **main switch**



2. Turn on **2 switches** in the rack (System + components)
3. Turn **the key** (laser I)



4. Start the PC and Zen software

Email: imaging@biomed.au.dk

Phone:

Nina: +45 93 52 18 88

Anna: +45 22 47 50 35

Shut Down LSM 800

1. Clean objectives and cover the microscope
2. Turn off Zen and the PC
3. Turn **the key** (laser 0)
4. Turn off **2 switches** in the rack (System + components)
5. Switch off the **main switch**
6. Note time of use and system status in the logbook

Email: imaging@biomed.au.dk

Phone:

Nina: +45 93 52 18 88

Anna: +45 22 47 50 35