


# Start Up LSM 900

- 1,2. Turn on 2 switches in the rack (System + components)
3. Turn the key to the position I (laser I)

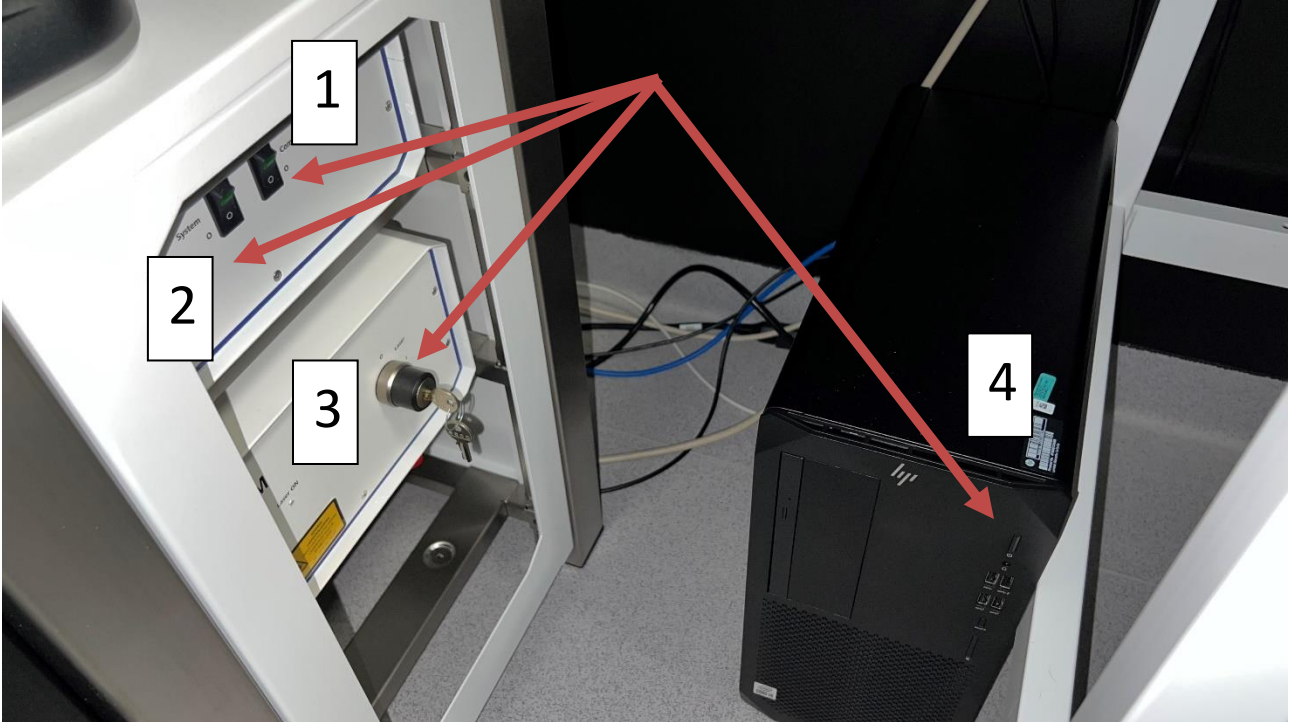

4. Start the PC and Zen software

# Shut Down LSM 900

1. Turn off Zen software
2. Transfer YOUR DATA from D: drive to O: drive.
2. Clean objectives and cover the microscope
2. Turn off the PC (4)
3. Turn the key to position 0 (laser 0) (3)
4. Turn off 2 switches (1,2) in the rack (System + components)
6. Note time of use and system status in the logbook

Email: [imaging@biomed.au.dk](mailto:imaging@biomed.au.dk)