

Start Up LSM 710

1. Turn on **3 switches** (Main switch, PC system, Components)



2. Turn on the **red switch** on the wall (Real Time PC)



3. Turn on Argon laser on by the **black switch** and turn **the key** (laser room)
(only if 458nm, 488nm and 514nm excitation is needed)



ON/OFF button Argon laser

ON/OFF Key Argon laser

4. Flip the **toggle switch** to "Run"



Toggle switch (up=run; down=standby)

Please don't touch "light control", but the correct position is just before the diode goes from green to red

5. Start the PC and Zen software
*NOTE!!! If Zen does not start up (activate boot status; restart Real time PC by **red switch** (point 2 above) and restart Zen*

Shut Down LSM 710

1. Flip **toggle switch** down



Toggle switch (up=run; down=standby)

2. Turn **key** Argon laser
3. Wait 10 min while lOaser cools down! Save data on your person drive while waiting.
4. Turn off **3 switches** (Main switch, PC system, Components)
Turn off the **red switch** on the wall (Real Time PC)
Turn off Argon laser on by the **black switch** and turn **the key**



ON/OFF button Argon laser

ON/OFF Key Argon laser

5. Shut down PC